

SIRCHIE®

FINGER PRINT LABORATORIES




Dazzle Fluorescent Magnetic Latent Print Powders

SIRCHIE's new line of Dazzle Fluorescent Magnetic Latent Print Powders offer a unique formulation for dusting latent prints. When dealing with confusing or multi-colored backgrounds, these fluorescent powders serve a dual purpose. The dark pigment of the Dazzle Powders is easy to discern under natural light. Depending on the Dazzle Powder used, they give off an intense fluorescent color of ORANGE, RED or YELLOW when excited.

They can be applied with all SIRCHIE magnetic powder applicators. View dusting area with an ultraviolet, blue or green forensic light source to see results.


DAZZLE ORANGE

The **FMP02** is naturally extreme dark orange, but fluoresces an intense orange when excited by UV light.

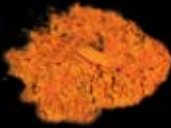

DAZZLE RED

The **FMP03** is naturally extreme dark red, but fluoresces an intense red.

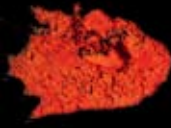

DAZZLE YELLOW

The **FMP04** is naturally extreme dark yellow, but fluoresces an intense yellow.



No. FMP03 Dazzle Red is used to develop latent prints on the multicolored surface of a magazine cover.

CAT. NO.	DESCRIPTION	PRICE
FMP02	Dazzle Orange, 1 oz. (30ml)	\$7.95
FMP03	Dazzle Red, 1 oz. (30ml)	\$7.95
FMP04	Dazzle Yellow, 1 oz. (30ml)	\$7.95

MAGNETIC POWDER APPLICATORS



No. 125L Standard Magnetic Powder Applicator



No. 125MD MEGAWAND™ Applicator



No. 125MBA Magnetic Burnishing Applicator...
its use with lifting material is shown below.



Standard Magnetic Applicator

The **125L** has the capacity to hold 25% more powder than ever before and weighs a mere 0.5 oz. (14.2g)! It is ideal for precision dusting of small to medium surface areas. The injection molded applicator features a narrow magnet to fit any size powder container. Process paper, finished and unfinished wood, nonferrous metals, plastic, glass, vinyl, etc.

MEGAWAND™ Magnetic Powder Applicator

The **125MD** has the capacity to hold 25% more powder than before. It features an expanded coverage area designed especially for dusting a large surface area. The wand holds approximately three times as much powder as the standard wand and features a machined Delrin® housing with spring-loaded ejection. Weight: 1.6 oz. (45.4g).

Magnetic Burnishing Applicator

The **125MBA** is capable of carrying more than three times the powder of standard wands and features an innovative burnishing head to improve the quality of crime scene latent lifts. Like its predecessor, the MEGAWAND™, the **125MBA** features a machined Delrin® housing with spring-loaded ejection. The difference is the built-in burnishing tool. Simply use the burnishing head to force out any air bubbles and wrinkles from the lifting material. Weight: 1.7 oz. (48.2g).



No. 125L

No. 125MD

No. 125MBA

CAT. NO.	DESCRIPTION	PRICE
125L	Standard Magnetic Powder Applicator	\$18.50
125MD	MEGAWAND™ Magnetic Powder Applicator	\$21.75
125MBA	Magnetic Burnishing Applicator	\$35.95